

# pH 7.00

## Calibration solution

**Works with any pH Probe**


**20ml each pouch**

**25 pouches per box**

**±0.02 pH at 25°C**

**Standardized against  
NIST-certified references**

**2 year shelf life at room temperature**

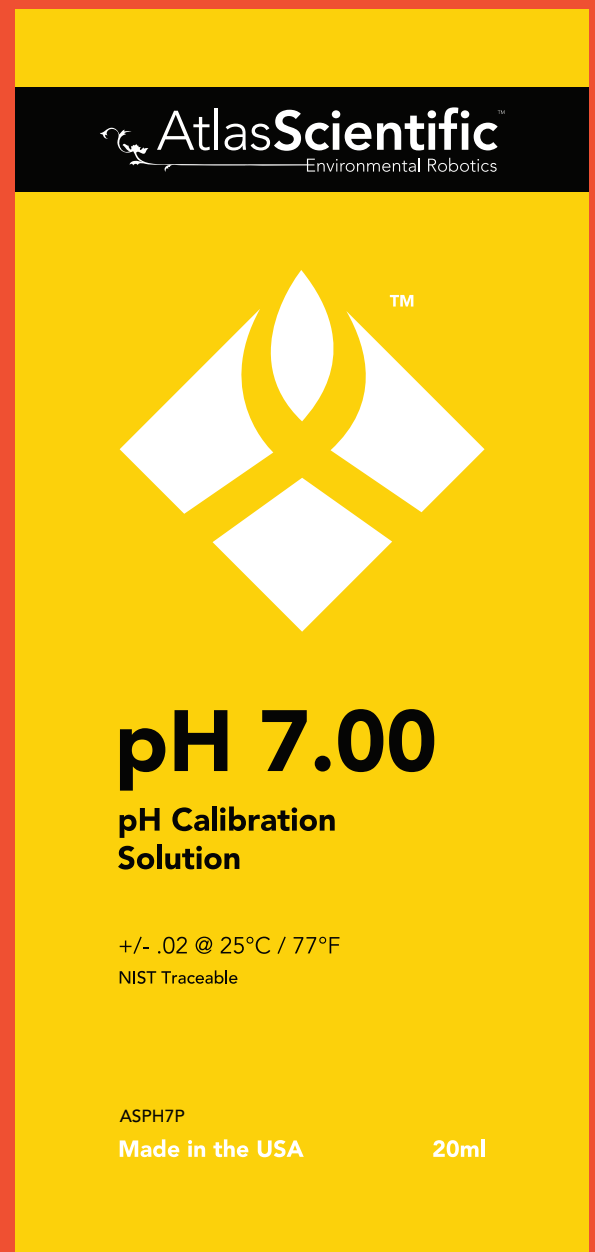


# Pre calibration

Remove the soaker bottle and rinse off the pH probe. Once the probe has been thoroughly rinsed, dry off the probe with a paper towel.



Remove the top section of the pouch and insert your pH probe. Let the probe sit in the calibration solution until readings stabilize.

This process should take about 1 – 2 minutes.

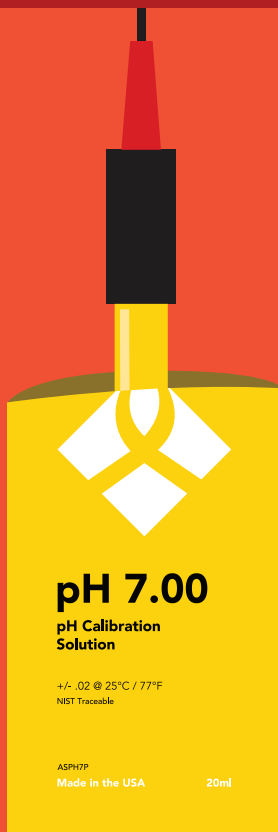

# 1 – 2 mins

Once readings have stabilized you are now ready to calibrate.

# Post calibration

Once the calibration is complete, dispose of the used solution.

